


	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

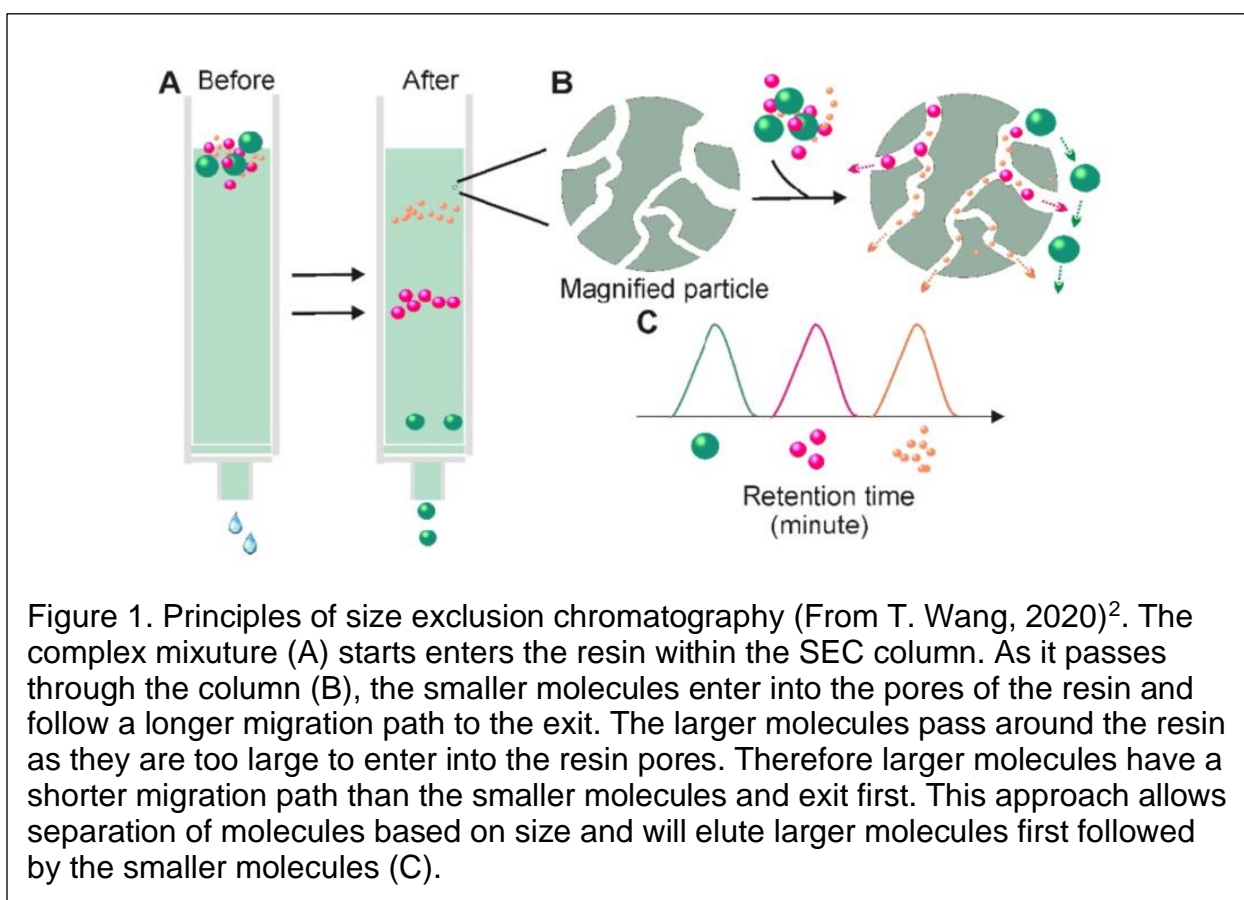
Background

Biological systems are complex. It is difficult to study the structure and function of individual molecular components of these systems or how they might interact with other components of the system. To overcome this challenge, scientists have developed a variety of methods to isolate, or purify, protein of a single type. These purified proteins are obtained as homogeneous aqueous solutions of the pure protein often stored in a compatible buffering solution to maintain protein solubility and stability. Methods of protein purification vary greatly and depend on the biochemical properties of the protein of interest. In this series of SOPs (“Protein Purification 1-3”), we will provide training and instruction on purifying protein through (1) ion exchange, (2) size exclusion chromatography, and (3) affinity purification.

The choice of which protein purification method to use depends on the biochemical properties of the protein of interest, as well as the desired degree of purity. The biochemical characteristics of a protein that can be used to allow for separation from other molecular components include its size, charge, stability (precipitation or solubility under various conditions), presence of any affinity tags, etc. When one decides to purify a protein, the above conditions are important to consider, but even knowing all of them, there can still be variability in approach depending on the protein. If the protein of interest is one that has been purified before, it is recommended to study past protocols proven effective and use or modify them as needed to suit your needs.



Our protein of interest is 1GFP, which is composed of the fusion protein His6-GFP-TEV. His6 is an affinity tag made up of six consecutive histidine residues. This affinity tag is useful for affinity purification using nickel (Ni^{2+}) chelated resin. GFP is our protein target of interest, a highly stable fluorescent beta-barreled protein that was first isolated from the jellyfish *Aequorea Victoria*.¹ The Tobacco Etch Virus (TEV) protease site is a small peptide sequence that is useful for recognition and cleavage by the TEV protease. If the TEV site was between the His tag and the protein of interest, this site would allow for controlled removal of the His6 affinity tag post expression or purification. This is an approach many use, but is not used here, nor is it needed for our purposes.

	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>



Together, this fusion protein product can be expressed, and purified by a number of different approaches, three of which are described in this series of SOPs.

In this protocol, we describe purification of 1GFP via size exclusion chromatography (SEC). SEC allows users to separate small molecules based on their size. To accomplish this, SEC columns are used in conjunction with the normal workflow of protein purification via HPLC. SEC columns contain a unique type of resin in which each bead of the resin contains small pores that small molecules can enter and pass through (Figure 1). This principle allows for separation of molecules based on size



	<p align="center">SKILLS CENTER</p> <p align="center">STANDARD OPERATING PROCEDURE</p>	<p align="center">A BIOFIZZ</p>  <p align="center">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

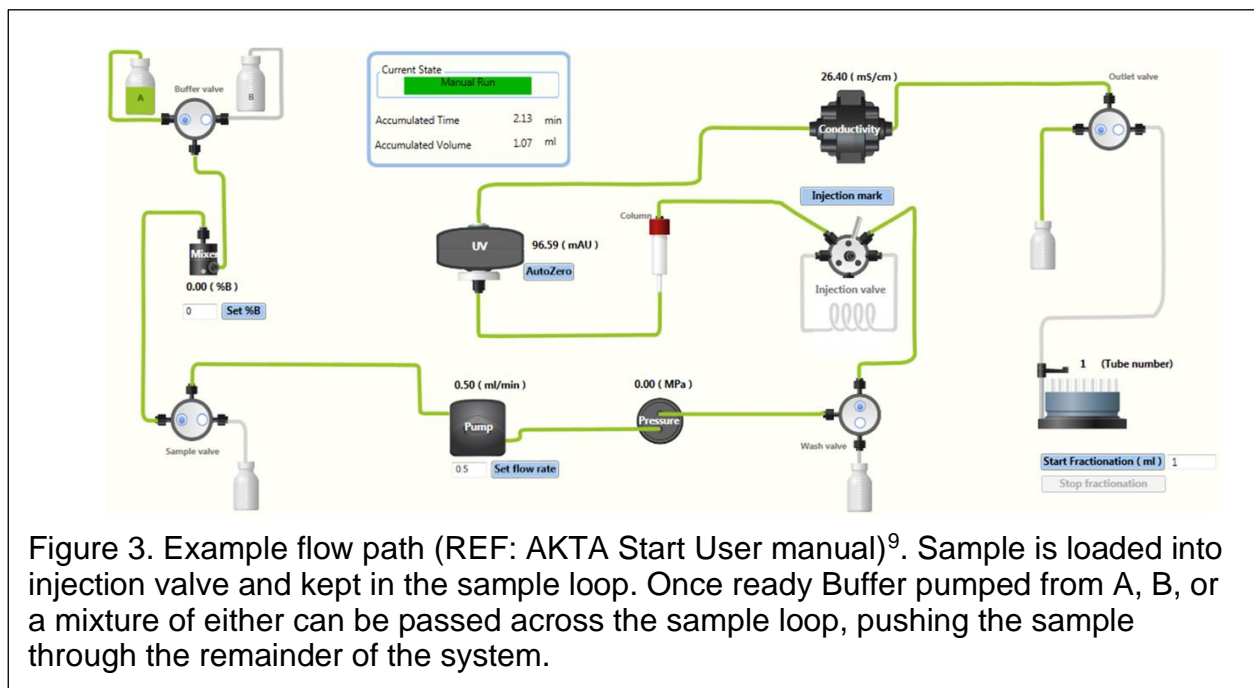
alone and does not select for differences in charge or affinity tags as the past methods have.

While this approach is simple and only requires one buffer, it is more likely to experience sample loss, compared to the previously described methods, as there is a longer path through resin required for the protein to travel. This method can also take longer to accomplish as protein purification using SEC columns often requires much slower flow speeds to maintain the structural integrity of the resin.



This approach can also be useful for exchanging buffers of your protein sample of interest. Loading the sample and using a different and new sample buffer can allow for collecting your sample at the end in the new buffer of interest. Though again, this comes at the risk of sample loss. The remainder of this SOP will instruct you in protein purification via SEC.



	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher



Each with an entry and exit valve, all of the protein purification columns described in this series of SOPs are connected to a machine that allows one to control the purification of the proteins across these ion exchange columns. These machines are known as **high-performance liquid chromatography (HPLC)** purification systems (Fig. 2). Aside from the HPLC purification columns, Figure 2B shows a diagram of the additional primary components of a HPLC system. These include the buffering solvent or solvents, a pump to control flow speed of the sample and buffer solutions, a small port before the HPLC column to allow for injection of the protein sample to be purified, a detector for monitoring separation of molecular components (usually in the form of a UV spectrophotometer), a fractionator (Turntable shown in Fig. 2 and 3) used to collect defined small volume aliquots of fractionated sample, and finally the waste container to collect the components not needed. The HPLC is connected to a computer that serves as the primary interface for operating the HPLC as well as collecting post column absorption data and the location of collected fractions across that data. As we will see

	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

more in later SOPs within this series, the HPLC serves as a highly versatile tool for controlled purification of liquid soluble proteins.

Given the quantity of components of HPLC protein purification and the information required to operate these types of systems, this skill can take time to master.. There are many small details to consider and questions you will generate as you familiarize yourself with the HPLC protein purification system. Please use your available resources (proctors, AKTA user manuals, etc) to ensure you are properly prepared before attempting to purify your protein of interest.

After one successfully purifies their protein across a particular chromatography column, it is common practice to run a series of samples of your collected fractions over SDS-PAGE to determine which samples are the are the purest and which you desire to collect, pool and store as your final product. After completion of the above, it is important to restore the HPLC system back to its default state with the proper wash cycles and filling of the lines with the proper storage buffer.

1. Purpose

The purpose of this SOP is to instruct students on the background and procedures necessary for protein purification via size exchange chromatography.

2. Scope

This procedure applies to qualified skills center users.



3. Responsibility

3.1. It is the responsibility of the user to understand and perform the procedure described in this document.

3.2. It is the responsibility of the user performing the procedure to fully document any deviations from the written procedure.

3.3. It is the responsibility of the user to become trained on the procedure.

4. Definitions

	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher



4.1 Chromatography – The process used to separate components of a soluble mixture. The mixture to be purified is composed of the mobile phase, which passes through the purification chromatography resin, and a stationary phase, which is captured or bound to the purification resin. Immobilized proteins of the stationary phase can be eluted using chemical components that stimulate the unbinding of protein from resin. Eluted protein is collected and stored as the purified final product.

5. Materials/Equipment

- 5.1 Ice bucket + ice
- 5.2 Harvested bacterial pellet of cells after expression of 1GFP.
 - 5.2.1 Obtained from the end of the protein expression SOP
- 5.3 Freshly prepared Lysis Buffer, Make lysis buffer if not available
 - 5.3.1 Liquid TritonX-100
 - 5.3.2 Powdered Sodium Phosphate Monobasic, Monohydrate (MW=137.99 Da)
 - 5.3.3 Powdered Sodium Phosphate Dibasic, Anhydrous (MW=141.96 Da)
- 5.4 Buffer A (See recipes below)
 - 5.4.1 Powdered Bis-Tris (MW=209.2418 Da)
 - 5.4.2 Powdered NaCl (MW=58.44 Da)
 - 5.4.3 500 ml graduated cylinder
 - 5.4.4 Stir bar (that fits base of graduated cylinder) + stir plate
 - 5.4.5 pH meter
 - 5.4.6 Concentrated HCL
 - 5.4.7 Glass filter sterilizer
 - 5.4.8 2x Clean autoclaved 500 ml buffer bottles

6. Recipes

- 6.1 500 ml Lysis Buffer: 10 mM NaCl, 1.0% TritonX-100, 50 mM Phosphate, pH 7.4
 - Measure 400 ml dH₂O in clean 500 ml graduate beaker
 - Add .292 g of NaCl
 - Add 5 mL of 100% TritonX-100
 - Add 3.55 g of Sodium Phosphate Dibasic, Anhydrous
 - Add 3.45 g of Sodium Phosphate Monobasic, Monohydrate
 - Bring up to total volume of 500 ml after dry reagents are dissolved
 - Adjust desired pH using concentrated HCl/NaOH

	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher

6.2 750 ml Buffer A: 20 mM bis-Tris buffer, 150 mM NaCl, pH 6.7.

Measure 600 ml dH₂O in clean 1000 ml graduated cylinder

Add 3.13 g powdered Bis-Tris

Add 6.57 g NaCl

Mix with stir bar on stir plate till dissolved

Starting pH will be ≥ 6.7. Measure pH and if needed, adjust to desired pH using concentrated HCL to final pH = 6.7.

Add dH₂O to final volume of 750 ml.

Filter sterilize.

REF: "How to Filter Buffers (Vacuum Filtration)" Training video → <https://www.youtube.com/watch?v=Haw14QOFxL8>

NOTE: Instructor mentions turning on "water." This is just the valve that provides vacuum suction from the tube connected to the filtering flask.

Add filter sterilized buffer to clean autoclaved 1000 ml flask.

Label with buffer, pH, date made, your name.

Can store at room temperature with other buffers.

7. Procedures (Adapted from Van Dyke M., 2021¹⁰)

7.1 Obtain bacterial pellet of harvested cells post induced expression of 1GFP.

7.1.1 Thaw slowly on ice and briefly with gloved hands to warm.



7.1.2 If you do not have a pellet of cells containing expressed protein, please stop and start at the protein expression SOP.

7.2 Measure approximate volume of thawed cells. Transfer all into 2 ml microcentrifuge tubes.

7.2.1 Save 100 µl aliquot of freeze/thawed sample (F/T). Label and store for later.

7.3 Resuspend the cell pellet in TX-100 Lysis buffer by pipetting or vortexing briefly until the suspension is homogenous:

7.3.1 Use 0.025 - 0.075 mL of Tx-100 Lysis buffer for every 1 UOD600 harvested. To calculate the UOD600, multiply the volume harvested by the OD600 reading.

	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher

-For example, a 5 mL culture harvested at OD600 1 gives 5 mL x 1.0 = 5 UOD600. In this example, 0.125 – 0.375 mL Lysis Reagent is required to lyse efficiently.

-If harvested cells are weighed, use 10 mL of lysis buffer per 1 gram of cells.

7.3.2 Incubate the resuspended cells at room temperature for 10 - 20 min with gentle shaking, gentle rotation, or swirling. Lysis is usually visible with a clearance of the suspension.

7.3.3 Save 150 µl aliquot of soluble sample. Label (“L/S”). Store on ice.

7.3.4 Dilute one of the 150 ul aliquots with 1.35 mL of Buffer A to be purified by IEX purification to make a total of 1.5 mL of sample in solution. Vortex Briefly.

7.3.5 Save 100 µl aliquot of lysis/soluble sample. Label (“L/S”). Store on ice.

7.4 Purify protein via size exclusion chromatography.

7.4.1 Prepare appropriate buffer. See recipe for preparation of buffer A above.

7.4.2 Be sure to familiarize yourself with the primary components of our lab’s HPLC (Fig. 2).

7.4.2 Familiarize yourself with the SEC column you will use → HisTrap excel, 1 ml (<https://cdn.cytivalifesciences.com/api/public/content/digi-11639-pdf>)



7.4.2.1 Document the following: Type of column, volume of column, max pressure limit, recommended pressure limit, max flow rate, recommended flow rate.

7.4.3 Make sure to always be aware of the flow path of buffers through the solution.

7.4.4 The column should already be connected to the HPLC and stored in EtOH. Check to ensure it is connected. Make note of the valves the column is connected to. We will tell the computer to direct the pathflow out these valves and across our column soon.

7.4.5 Prepare HPLC, column, and fractionator

7.4.5.1 Pump wash both pumps A & B.

	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher

-Make sure to open the wash valve and plug the top of the mixer avoid washing the column with high flowrate buffer

7.4.5.2 Set high pressure alarm, set flow speed, and select the pump that will pump water through the system.

7.4.5.3 Set UV detector to 280 nm (for protein) and 255 nm (for detecting potential nucleic acid contaminants).

7.4.5.3 Flow 1.5 column volumes (CVs) of water through your column. This will displace the EtOH it was stored in.

7.4.5.4 Remove tubing lead from water (weighted ends of HPLC tubing currently in water flasks; Should be labelled "A" for the pump A). No need to rinse leads as they were previously in water. Place "A" tubing lead into buffer A.

7.4.5.5 Pump wash with buffer solutions. This prepares your pumps with the buffer that will soon pass over your column.

7.4.5.6 Wash column as in previous step with buffer A. Make sure that the baseline UV reading is flat after low salt buffer wash. This is your baseline. Presence of your protein eluting from the column will be observed with noticeably large 280 nm peak.

7.4.5.7 Make sure fractionator is in place.

7.4.5.8 Load fractionator with appropriate collection tubes.

If collecting up to 1 ml samples, you can fill fractionator with full sized collection tubes with 1.5 ml capless epi tubes sitting within each large tube.

7.4.5.9 Fill rack with 65 collection tubes.



7.4.6.10 Lastly, place the waste output tubing into a large clean empty flask to collect waste.

7.4.6 Prepare protein sample

7.4.6.1 Visually inspect protein sample. If you observe large quantities of precipitate, centrifuge briefly (high speed, 5 min, 4°C), collect supernatant and transfer to new tube.

7.4.6.2 Filter protein containing sample through Millex-HP 0.45 µm pore sized syringe filter into a clean receptacle.

7.4.6.3 Take note of the location, and volume of the sample loop (See figure 4), the looping of tubing sized to hold your protein sample prior to its

	<p align="center">SKILLS CENTER</p> <p align="center">STANDARD OPERATING PROCEDURE</p>	<p align="center">A BIOFIZZ</p>  <p align="center">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

injection onto the column. Make sure the volume of your sample loop is greater than, or equal to the total volume of your sample. If the sample loop is smaller than the total volume of your sample, exchange it for an appropriately sized sample loop.


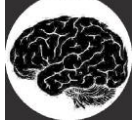
7.4.6.3 Select a glass syringe with a volume large enough to hold your entire sample.

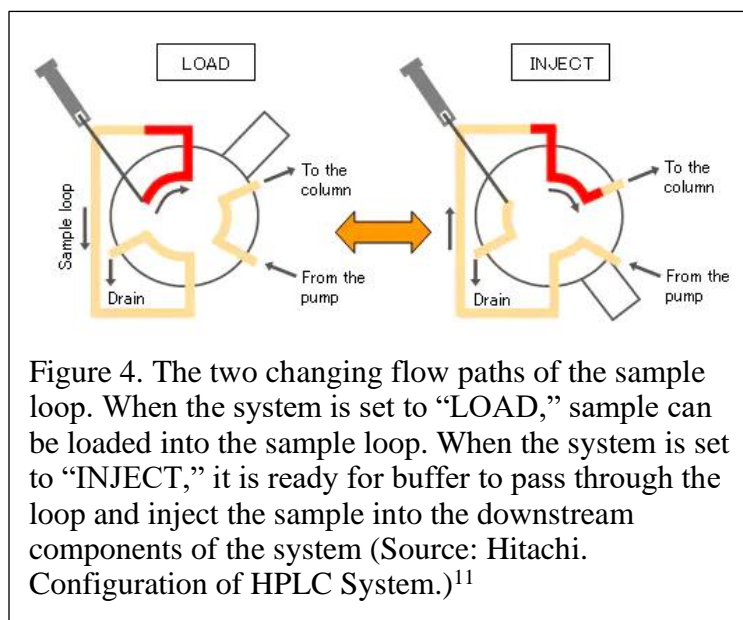
7.4.6.4 Carefully draw up the filtered sample into a glass syringe. Carefully displace any air bubbles.

NOTE: Introducing air bubbles into the HPLC and connected columns has the potential to disrupt the viability of the column and proper functioning of the HPLC. Please exercise caution to prevent this. Contact a proctor if you have any concerns regarding this.



7.4.7 Load protein sample onto sample loop (Read all prior to execution)

7.4.6.5 Before injecting, place syringe tip into injection port, resting on the injection port syringe holder.

	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>



- 7.4.6.6 The injection valve should be preset to “LOAD.” Make sure to check this setting in the manual injector settings. THEN you are free to load your sample into the super loop. Leave syringe connected to injection valve and resting on syringe holder.
- 7.4.8 Double check everything prior to run.
- 7.4.8.1 Ensure proper preparation of HPLC and buffers, full fractionator to collect samples, and empty flask to collect waste.
- 7.4.9 Run pre-prepared **SEC protein purification run**
- 7.4.9.1 Select pre-prepared purification protocol.
- 7.4.9.2 Make sure high pressure alarms match those you have been using.
- 7.4.9.3 Select “Run.”
- 7.4.9.4 Monitor the beginning stages of HPLC protein purification. Make sure the samples begin to be dispensed into the fractionator collection tubes, and that no pressure alarms go off.
- 7.4.10 AFTER RUN IS COMPLETE, observe the UV absorption data for your run. You should have a strong peak at 280 nm indicative of your purified protein.

	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

You will likely have other smaller peaks that come from non-target protein being passed through the UV detector. Ensure that peaks from other absorbances do not correlate with your protein's 280 nm peak. If there are, then you might still have contamination of other molecules in your purified protein sample. Other protein purification steps can be taken to further purify your protein if needed.

7.4.10.1 Cover your fraction collector loosely with seran wrap.

7.4.10.1 Print a copy of the UV absorbance data.

7.4.10.2 Select a sampling of fractions that capture the beginning, middle, and end of prominent peaks, as well as a few others within each peak. Mark these on your printed UV data sheet.

7.4.10.3 Gather enough 0.5 ml epi tubes to capture all peaks you would like to test for protein.

7.4.10.4 Fill each with 2 μ l 6X SDS-PAGE running buffer (dark blue).

Label these tubes with the corresponding fraction numbers you will test for protein.

7.4.10.5 Add 10 μ l of each sample you've selected into the corresponding tube containing SDS-PAGE running buffer. Gently mix via pipetting.

7.4.10.6 Recover fractionator containing samples. Temporarily store in 4°C fridge. Can be stored overnight.

7.4.11 Along with an appropriate SDS-PAGE ladder, and all labelled pre-purification samples, run all samples over SDS-PAGE and stain with Coomassie. Consult SDS-PAGE SOP for details if needed.

7.4.11.1 After gel is de-stained, observe bands that you believe to be corresponding to the molecular weight of your protein of interest.



Document notes.

7.4.12 Pool all samples corresponding to the most-pure fractions of your protein of interest. Dispose of the rest.

7.4.12.1 Measure concentration of purified protein sample via UV VIS spectrophotometer.

Consult proctors for help if needed.

7.4.12.2 Measure your purified protein's UV Vis *Spectra* (Y-axis: Absorption, X-axis: Wavelength). Compare spectra to your observations about the observed protein's color. Why do you think the sample is green?

	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher

7.4.12.2 Label sample of pure protein with your name, date, the protein, the concentration of protein, and the contents and pH of the buffer the protein is in.

7.4.16 In many applications, proteins can be concentrated for their specific use if needed after this final step.

7.4.13 Clean up

7.4.13.1 Wash HPLC

7.4.13.1.1 Set pressure alarms for all below. Run 2+ CVs of buffer across column.

7.4.13.1.2 Remove weighted buffer tubing leads from buffers. Rinse ends of leads and tubing with DI from a squirt bottle over an unused beaker.

7.4.13.1.2 Return leads to DI water flask.

7.4.13.1.3 Run water through the column.

7.4.13.1.4 Lower flow speed (as EtOH is more viscous than water) and run EtOH through system and through the column. The column can be stored in EtOH.

7.4.13.2 Rinse Super loop

7.4.13.2.1 Rinse glass syringe with clean DI water three times

7.4.13.2.2 Using this glass syringe or other clean syringe, rinse super loop with three or more super loop volumes with DI water.

7.4.13.3 Empty fraction collector and dispose of waste

7.6.13.3.1 0.5 ml capless epi tubes used to collect fractions that now do not contain your protein of interest can be disposed of.



7.6.13.3.2 Any larger fraction collection tubes can be rinsed a few times with DI water.

7.4.13.4 Dispose of any unused buffers.

7.4.13.5 Clean up any remaining used space in the lab.

8. Troubleshooting

8.1 There are many places for complications in purifying protein via a HPLC machine. Please read through notes carefully prior to executing protocol. Contact available proctor for help as needed.



	SKILLS CENTER STANDARD OPERATING PROCEDURE	A BIOFIZZ  PRODUCTON
Protein Purification (SEC) Module Hours: 5	Effective Date: 11/3/2023 PRQs: Protein Expression	Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher

9. References

1. Zimmer M. *GFP: from jellyfish to the Nobel prize and beyond*. Chem Soc Rev. 2009 Oct;38(10):2823-32. doi: 10.1039/b904023d. Epub 2009 Jun 15. PMID: 19771329.
2. Yang, Dongbin et al. (2020). Progress, opportunity, and perspective on exosome isolation - Efforts for efficient exosome-based theranostics. *Theranostics*. 10. 3684-3707. 10.7150/thno.41580.
3. Aryal S., 2022. *HPLC- Definition, Principle, Parts, Types, Uses, Diagram*. Microbe Notes – Instrumentation. <https://microbenotes.com/high-performance-liquid-chromatography-hplc/>
4. Laboratory Notes. Preparation of lysozyme stock solution (10 mg/ml). <https://www.laboratorynotes.com/preparation-of-lysozyme-stock-solution-10-mg-ml/>
5. AKTA Start User manual. https://bio.pnpi.nrcki.ru/wp-content/uploads/2020/01/GE_AKTA_start.pdf
6. M. Van Dyke. *Expression and Purification of Thermostable Proteins Expressed in E. coli*. 2021. dx.doi.org/10.17504/protocols.io.bq4jmyun
7. Hitachi. Principle and System Configuration of HPLC. <https://www.hitachi-hightech.com/global/products/science/tech/ana/lc/basic/course3.html>

10. MODULE METHODS TASK

- 10.1 Starting at the post-expression bacterial pellet, purify 1GFP via size exclusion chromatography.
- 10.2 Describe how, specifically, SEC can be used to purify our protein of interest.
- 10.3 What are the variables we must consider when choosing this as our protein purification method?
- 10.4 Display your results of the UV absorbance (A280) over time/volume from your purification run. Clearly label which peaks you believe to be corresponding to your protein of interest. Label which fractions you used to test for the presence of your protein.
- 10.5 Display your results from your SDS-PAGE gel run from all collected samples. Identify which bands correspond to your protein of interest, and why you believe those bands are your desired protein.

	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

- 10.6 Attach a photo of your final purified protein product containing the tube labelled with your name, date, protein, protein concentration, and buffer information.
- 10.7 Do you have any evidence that your protein sample is not completely pure? If so, what evidence does this consist of? What steps might you take next to further purify your protein of interest? What is a drawback of subjecting your protein to an entirely separate purification step or steps?
- 10.8 Please provide us with any feedback you wish to offer: What you liked about this training module and what you feel would be helpful to change.

11. 1GFP Information

11.1.1 Plasmid vector: pET His6 GFP TEV LIC cloning vector

11.1.2 Gene of interest: His6 GFP TEV (1GFP)



11.1.3 Gene length: 801 bp

11.1.4 Gene sequence:

5'

**ATGGGTTCTTCTCACCATCACCATCACCATGGTTCTTCTGTGAGCAAGGGCGAG
GAGCTGTTACCGGGGTGGTGCCCATCCTGGTCGAGCTGGACGGCGACGTAA
ACGGCCACAAGTTCAGCGTGCGCGGCGAGGGCGAGGGCGATGCCACCAACG
GCAAGCTGACCCTGAAGTTCATCTGCACCACCGGCAAGCTGCCCGTGCCCTG
GCCACCCCTCGTGACCACCCTGACCTACGGCGTGCAGTGCTTCAGCCGCTACC
CCGACCACATGAAGCAGCAGACTTCTTCAAGTCCGCCATGCCCGAAGGCTAC
GTCCAGGAGCGCACCATCTCCTTCAAGGACGACGGCACCTACAAGACCCGCG
CCGAGGTGAAGTTCGAGGGCGACACCCTGGTGAACCGCATCGAGCTGAAGGG
CATCGACTTCAAGGAGGACGGCAACATCCTGGGGCACAAGCTGGAGTACAAC
TCAACAGCCACAACGTCTATATCACGGCCGACAAGCAGAAGAACGGCATCAAG
GCGAACTTCAAGATCCGCCACAACGTCGAGGACGGCAGCGTGCAGCTCGCCG
ACCACTACCAGCAGAACACCCCATCGGCGACGGCCCCGTGCTGCTGCCCGA
CAACCACTACCTGAGCACCCAGTCCAAGCTGAGCAAAGACCCCAACGAGAAGC
GCGATCACATGGTCCTGCTGGAGTTCGTGACCGCCCGGGATCACTCTCGG
CATGGACGAGCTGTACAAGGGGATCGAGGAAAACCTGTACTTCCAATCCAATA
TTGGAAGTGGATAA**

-3'

	<p style="text-align: center;">SKILLS CENTER</p> <p style="text-align: center;">STANDARD OPERATING PROCEDURE</p>	<p style="text-align: center;">A BIOFIZZ</p>  <p style="text-align: center;">PRODUCTON</p>
<p>Protein Purification (SEC)</p> <p>Module Hours: 5</p>	<p>Effective Date: 11/3/2023</p> <p>PRQs: Protein Expression</p>	<p>Revision 1.0 Author: Zachary Hazlett, Oscar Cordova Checked by Editor: Carson Butcher</p>

11.1.5 Peptide sequence:

N-term

MGSSHHHHHHGSSSVSKGEELFTGVVPILVELDGDVNGHKFSVRGEGEGDATNG
KLTLKFICTTGKLPVPWPTLVTTLTLYGVQCFSRYPDHMKQHDFFKSAMPEGYVQ
ERTISFKDDGTYKTRAEVKFEGDTLVNRIELKGIDFKEDGNILGHKLEYNFNHSHNV
YITADKQKNGIKANFKIRHNVEDGSVQLADHYQQNTPIGDGPVLLPDPNHYLSTQS
KLSKDPNEKRDHMLLEFVTAAGITLGMDELYKGIENLYFQSNIGSG

C-term

11.1.6 Protein Parameters (ProtParam)

Expasy: <https://web.expasy.org/cgi-bin/protparam/protparam>

Number of amino acids: 266

Molecular weight: 29.752 kDa

Theoretical Isoelectric point: pI = 5.71